CONGEN

**SureFood[®] ANIMAL ID Pork
SENS PLUS**

Art. No. S6017
100 rxn

User Manual

January 2019

 Inhalt /  Content

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1 General Information

1.1 Description

The test detects pork DNA (*Sus scrofa*). Each reaction contains an internal amplification control. The real-time PCR assay can be used with established real-time PCR instruments, equipped for detection of two fluorescence emissions at 510 nm and 580 nm (FAM and VIC/HEX) at the same time. The technical validation of instruments was performed on Agilent Mx3005P, Agilent AriaDx, BioRad CFX 96, Roche LightCycler® 480 II, Roche cobas z 480 Analyzer, Roche LightCycler® 2.0², Applied Biosystems 7500, LTF MyGo Pro, R-BIOPHARM RIDA®CYCLER and Qiagen RotorGene Q.

1.2 Limit of Detection

The SureFood® ANIMAL ID Pork SENS PLUS real-time PCR is developed for the detection of pork-DNA in muscle meat mixtures at a relative amount of ≤ 0.0001 %. Due to the high sensitivity the method is suitable for the analysis of highly processed products like e.g. gelatin, sweets or meat bone meal. The assay limit of detection depends on sample matrix, processing grade, DNA preparation and DNA content.

1.3 DNA-preparation

For DNA-preparation the use of SureFood® PREP Basic and for highly processed food and feed the use of SureFood® PREP Advanced is recommended.

1.4 Kit components and storage

Kit Code	Reagent	Amount	Lid Color
1	Reaction Mix	2 x 1100 µl	Yellow
2	Taq Polymerase	1 x 11 µl	Red
3	Positive Control	1 x 200 µl	Light Blue

Store all reagents at -20°C and protected from light.

1.5 Additionally required equipment and materials

- DNA-Extraction kit (e.g. SureFood® PREP Basic or SureFood® PREP Advanced)
- real-time PCR instrument with two detection channels (510 nm, 580 nm)
- real-time PCR consumable (plates, tubes, capillaries, foils, caps) pipettes with filter tips
- unpowdered disposable gloves
- Vortex mixer
- micro centrifuge with a rotor for the reaction tubes

²For the use of the Roche LightCycler® 2.0 a Color Compensation is necessary. The SureCC Color Compensation Kit II (Art. No. F4010) must be used for the color compensation of such devices.

1.6 Setup

	BlockCycler/ RIDA®CYCLER/ LightCycler® 480 II	Rotorgene
Initial Denaturation (HOLD)	5 min, 95°C	1 min, 95°C
Cycles	45	45
Denaturation	15 sec, 95°C	10 sec, 95°C
Annealing/Extension (CYCLE)	30 sec, 60°C	15 sec, 60°C
Temperature Transition Rate/ Ramp Rate	Maximum	Maximum
Fluorescence Detection Setup (exemplary)	Detection: End of Extension Phase <u>Detection System pork:</u> Various devices FAM-channel, Quencher: BHQ Rotor-Gene Q Green LightCycler® 2.0 530-none LC480 I 483 nm, 533 nm LC480 II 465 nm, 510 nm <u>Internal Amplification Control:</u> Various devices VIC/HEX-Channel, Quencher: BHQ Rotor-Gene Q Yellow LightCycler® 2.0 560-none LC480 I 523 nm (excitation) - 568 nm (emission) LC480 II 533 nm (excitation) - 580 nm (emission) Passive Reference: None	
Detailed information on the setup of several real-time PCR devices is available at the CONGEN homepage: http://www.congen.de/en/company/downloads		

2 Qualitative Analysis

2.1 Protocol

2.1.1 Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions). Recommended control reactions: negative control, positive control and extraction control. The master-mix includes an internal amplification control (inhibition control) for each reaction.

It is also recommended to prepare the master-mix with 10 % additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix by vortexing and centrifuge before opening and use. The tube of the Taq Polymerase should be kept at -20°C and not be mixed by vortexing.

Example for the calculation and preparation of 10 reactions:

Components for master-mix	Amount per reaction	10 reactions (with 10% excess)
Reaction Mix	19.9 µl	218.9 µl
Taq Polymerase	0.1 µl	1.1 µl
Total volume	20 µl	220 µl

Mix each master-mix well and centrifuge shortly before use.

January 2019

2.1.2 Preparation of the real-time PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells or capillaries.
- Close the tube of the negative control (the negative control is ready for PCR without any addition).
- Pipette 5 µl of sample DNA into the designated tubes/wells or capillaries.
- Pipette 5 µl of the Positive Control into the designated tubes/wells or capillaries.
- Centrifuge all tubes/wells or capillaries shortly at low speed.
- Place tubes/wells or capillaries into the real-time PCR instrument and start the run according to the setup.

2.2 Interpretation of results

The evaluation has to be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer.

The results of the control reactions have to be correct.

Pork DNA is detected in the FAM-channel. The internal amplification control will be detected in the VIC/HEX-channel.

A sample is stated **positive** for pork, if the sample DNA shows amplification in the FAM-channel. A sample is stated **negative** for pork, if the sample DNA shows no amplification in the FAM-channel and the internal signal (VIC-channel) of the sample is **positive**.

In the case of a negative result in all channels including the VIC-channel the sample contains PCR inhibiting substances. Under these circumstances an evaluation of the sample is not possible. DNA isolation and purification for the sample need to be improved.

3 Further Information**3.1 Product Information**

- Validation Report

3.2 Technical Support

For further questions please contact your distributor or send an e-mail to sales@r-biopharm.de.

3.3 Distribution and ordering

R-Biopharm AG
An der neuen Bergstrasse 17,
64297 Darmstadt, Germany
Phone: +49 (0) 61 51 - 81 02-0
Fax: +49 (0) 61 51 - 81 02-20
E-Mail: orders@r-biopharm.de
www.r-biopharm.com