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RIDA[®] Extraction Solution (colorless)

1. Intended Use

1.1. Use with RIDASCREEN[®]FAST Casein (R4612)

The RIDA[®] Extraction Solution (colorless) (R7098) can be used alternatively to RIDA[®] Extraction solution / Extraktions-Lösung (R7099) for the sample preparation of sausages and bakery goods in the RIDASCREEN[®]FAST Casein (R4612). R7099 is stained to simplify pipetting.

The sample preparation is described in the product information or application of the test.

1.2. With Gliadin test systems

The RIDA[®] Extraction Solution (colorless) (R7098) can be used for the sample preparation of raw materials, heated and processed food for the following qualitative tests:

- RIDA[®]QUICK Gliadin (R7003)
- RIDA[®]QUICK Gliadin (single packaged) (R7004)

It is recommended to use the Cocktail (patented) (R7006/R7016) for quantitative analysis with RIDASCREEN[®] Gliadin (R7001) or RIDASCREEN[®]FAST Gliadin (R7002).

The RIDA[®] Extraction Solution (colorless) (R7098) should only be used after comparison with Cocktail (patented) (R7006/R7016), because extraction with RIDA[®] Extraction Solution (colorless) (R7098) may result in reduced recoveries in some samples. An application can be requested at R-Biopharm.

2. Content

105 ml

3. Precautions and Handling

The RIDA[®] Extraction Solution (colorless) is harmful to health. It contains mercaptoethanol, therefore working under a chemical hood is recommended and skin contact should be avoided.

The solution should only be used by trained laboratory employees.

For hazard notes on the contained substances please refer to the appropriate material safety data sheets (MSDS) for this product, available online at www.r-biopharm.com.

4. Storage and disposal

Store at room temperature (20 - 25 °C / 68 - 77 °F).

Do not allow the product to reach sewage system and dispose of contents in accordance with national regulations.

The data corresponds to our present state of technology and provides information on our products and their uses. R-Biopharm makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. Defective products will be replaced. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. R-Biopharm shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.

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are registered trademarks of R-Biopharm AG