Compact Dry Swab

REF **ZCS1002953**

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1. Intended use

The Compact Dry Swab is a swab unit consistent of a sterile swab in a tube with 1 ml sterile PBS peptone water buffer for sampling, transport and subsequent quantitative detection of present microorganisms in combination with Compact Dry culture medium plates.

2. General information

The Compact Dry Swab is designed for sampling on surfaces in the production environment. The buffer contained in the tube is used as liquid stabilization and transport medium. The Compact Dry Swab enables hygiene control of surfaces using Compact Dry culture medium plates.

3. Test principle

The surface sample will be taken with the swab stick. The sample will be homogeneously mixed in the buffer medium by vigorously shaking. Subsequently the liquid has to be applied completely on a Compact dry plate using the outlet in the bottom of the tube. The inoculated culture plates have to be incubated according to their specification (see individual Compact Dry instructions for use) and evaluated by counting the number of specific colonies.

4. Reagents provided

In each package bag are 40 Compact Dry Swabs included.

Component	Cap color	Format	Volume
Swab unit	Orange	Ready to use	40 Pieces

Composition per swab unit:

- 1 Transport tube with outlet screw-cap
- 1 Swab stick attached to the orange colored screw-cap
- 1 ml PBS peptone water buffer (peptone 1.0 g/l, sodium chloride 5.0 g/l, potassium di-hydrogen phosphate 1.5 g/l, di-sodium phosphate (water free) 3.57 g/l, (9.0 g/l dodecahydrate), pH 7.0 ± 0.2)

The buffered peptone water (BPW, ISO 6887-1:2017) is prepared in accordance with the usage and is aseptically packaged.

The complete product is sterilized by gamma irradiation after packaging.

5. Required reagents / equipment (not provided)

- Incubator (30 - 35 °C / 86 - 95 °F): e.g. CULTURA® Mini-Incubator (Art. No. ZC7140651)
- Compact Dry plates (Art. No. HS7311 – HS9802)
- Optional: Frames for definition of 100 cm² (Art. No. ZFR1600000)

6. Warnings and precautions for the users

This test should be carried out only by trained laboratory employees. The instruction for use must be strictly followed.

The Compact Dry Swab is only suitable for in-vitro-use.

Please refer to the component safety information in the material safety data sheets (SDS) for this product, available online at www.r-biopharm.com.

Ensure the proper and responsible disposal of all reagents and materials after their use. For disposal, please adhere to national regulations.

7. Storage instructions

Store the Compact Dry Swab at room temperature (5 - 30 °C / 41 - 86 °F).

Do not use the test kit after the expiration date (see test kit label).

8. Indication of instability or deterioration of reagents

Tubes with stained or turbid liquid should be discarded.

9. Test procedure

1. Write the sample ID on the buffer tube using a water resistant pen.
2. Unscrew the orange swab stick from the tube.
3. Wipe the desired sampling surface (20 - 100 cm²; frames for definition of 100 cm², Art. No. ZFR1600000, available at R-Biopharm) with the swab and replace the stick back into the tube and screw it in tightly.
4. Homogenise the sample by shaking the tube.
5. Open outlet screw-cap and apply the whole solution content (1 ml) of the tube to a Compact Dry plate by pressing on the smooth middle part of the tube.
6. If necessary, prepare further series of 10 times sample dilution and inoculate 1 ml each in the middle of the Compact Dry plate.

7. Incubate the medium plates under specified conditions (see individual Compact Dry instructions under “Incubation time / Incubation temperature”).
8. Count the specific colonies which are grown.

Detailed instructions for the use of Compact Dry are included in the respective test kits.

Remark:

- Please do not touch the swab.
- Since the Compact Dry plates detects up to 3×10^2 colonies (cfu) per plate, original solution may be inoculated direct on Compact Dry plate without any dilution in case the expected number of bacteria is less than 3×10^2 cfu.
- The Compact Dry Swab and Compact Dry Plates are manufactured in an ISO 9001 certified facility.

An overview of available Compact Dry products, package sizes and article numbers is included in the Compact Dry product brochure and R-Biopharm product catalogue (available on request).

Further product information, please contact your local distributor or R-Biopharm at this address: sales@r-biopharm.de.

Version overview

Version number	Chapter and title
2016-02-08	Release version
2021-01-14	4. Reagents provided, 7. Storage instructions, 9. Test procedure + Remark

Explanation of symbols

- General symbols:

	Follow the instructions for use
	Batch number
	Expiry date (YYYY-MM)
	Storage temperature
	Article number
	Number of test determinations
	Manufacturing date (YYYY-MM)
	Manufacturer + address

Disclaimer

The user assumes all risk in using R-Biopharm AG's products and services.

R-Biopharm AG will warrant that its products and services meet all quality control standards set by R-Biopharm AG, and R-Biopharm AG will, at its option, replace or repair any components, product or repeat services which prove to be defective in workmanship or material within product specific warranty periods or expiration dates and which our examination shall disclose to our satisfaction to be defective as such.

This warranty is expressly in lieu of all other warranties, expressed or implied, as to quality, description, fitness for any particular purpose, merchantability, productiveness, or any other matter. R-Biopharm AG shall be in no way responsible for the proper use of its products and hereby disclaims all other remedies, warranties, guarantees or liabilities, expressed or implied, arising by law or otherwise, and it shall have no liability for any lost profits or damage, direct, indirect or otherwise, to person or property, in connection with the use of any of its products or services.

This warranty shall not be extended, altered or varied except by a written instrument signed by an authorized representative of R-Biopharm AG.

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are registered trademarks of R-Biopharm AG.