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# QuickGEN PCR Kit

## Alicyclobacillus Differentiation

### 1. Intended use

Qualitative detection and differentiation of *Alicyclobacillus* spp., *Alicyclobacillus acidocaldarius* and *Alicyclobacillus acidoterrestris* .

### 2. Test principle

The (TaqMan<sup>®</sup>) real-time PCR is based on hot-start-PCR and sequence-specific dual labelled probes (FAM/DQ; CY5/DQ; HEX/DQ) which, when accurately hybridised, emit a measurable fluorescent signal of a defined wavelength in the extension phase.

The increase of signal is continuously measured in a real-time PCR detection instrument. To avoid false negative PCR-results an Inhibition control (ROX/DQ) is amplified together in one reaction vessel with the specific sequence. The system contains dUTP. Optional: the use of Uracil-N-Glycosylase will eliminate any contamination with Uracil containing amplicons from former PCRs (the enzyme is not part of this kit).

### 3. Kit contents

The kit contains sufficient reagents for 48 PCR reactions:

1 x Premix	white cap
48 x Dye Strips (freeze-dried, incl. IC-DNA)	tube strips
6 x Cap Strips	
1 x ddH <sub>2</sub> O	colourless cap
1 x Control-DNA (freeze-dried)	yellow cap

## 5. Storage conditions

**The Control-DNA is freeze-dried and has to be solved in ddH<sub>2</sub>O prior to use (see 7.1).**

Do **not** freeze the lyophilized Dye Strips and lyophilized Control-DNA.

The PCR reagents should be stored at 2 – 8 °C (35 – 46 °F).

Keep Premix for storage at - 20 °C (- 4 °F) after opening. Avoid loss of sensitivity by repeating freezing and thawing more than 3 times. For irregular use aliquot the Premix.

The Dye Strips contain the fluorescent labelled probes and should be handled light protected. All reagents are stable for 12 months, if they are stored correctly.

## 5. Materials required, but not provided

### 5.1. Instruments

Real-time PCR machine for high profile PCR tubes

Standard benchtop mini-centrifuge (1.5 – 2.0 mL vials)

Centrifuge for strips

Pipettes

“Vortex”

### 5.2. Reagents and plastic ware

sterile ddH<sub>2</sub>O

sterile filter tips

**optional:** Uracil N-Glycosylase (0.01 U/μL added to the PCR reaction)

## 6. PCR

### 6.1. PCR-Setup

*When using the kit for the first time, the freeze-dried Control-DNA has to be shortly centrifuged and carefully resolved:*

- add 55  $\mu\text{L}$  sterile ddH<sub>2</sub>O to the freeze-dried Control-DNA
- after 15 minutes mix well

**Before every use thoroughly mix all PCR components and centrifuge briefly.**

Remove the film from the needed tube strips and pipette the PCR-Components. After pipetting close the tube strips with the provided cap strips.

PCR-reaction per sample:

PCR-Components	volume ( $\mu\text{L}$ )
Premix	15.0
Sample-DNA	5.0*
Total volume	<b>20.0</b>

\* if using the Simplex easy DNA kit: add 2.5  $\mu\text{L}$  DNA and 2.5  $\mu\text{L}$  sterile ddH<sub>2</sub>O

1. Pipette 15  $\mu\text{L}$  of the Premix into each PCR-tube, making sure that, prior to the first filling, the tip has been moistened.
2. Add 5  $\mu\text{L}$  sample DNA, add 5  $\mu\text{L}$  of the Control-DNA for the positive control reaction, add 5  $\mu\text{L}$  of the extraction control and 5  $\mu\text{L}$  of ddH<sub>2</sub>O for the negative control\* reaction. Use a fresh tip with each DNA filling.
3. Close the tubes immediately and centrifuge them shortly.
4. Place the tubes in the PCR-machine and start run.

**Very important: \* Please fill up the negative control with 5  $\mu\text{L}$  ddH<sub>2</sub>O to avoid unspecific amplification.**

**Work swiftly to avoid warming up and keep away from light.**

## 6.2 PCR-Program

The Premix contains no ROX. This must be considered according to the settings of the real-time machine. **ABI 7500:** Under “Assign Targets and Samples” in ‘Select the dye to use as the passive reference’ choose „none“.

**Mx3500P:** Under Instruments in ‘Filter Set Gain Settings’ change the Filter set gain multiplier to CY5 x1, ROX x1, HEX-JOE x4 und FAM x8.

For the use of UNG the thermal cycler program has to be changed according to manufacturers instructions.

Step	Time	Temp.	
Initial denaturation	15 min	95 °C	
Cycling Denaturation	15 sec	95 °C	Cycle 40 x
Cycling Annealing/ Elongation	25 sec	60 °C	

## 7. Evaluation

The evaluation has to be made according to the data analysis programme recommended by the real-time instrument manufacturer.

**Alicyclobacillus spp.-DNA:** HEX-channel

***Alicyclobacillus acidoterrestris*-DNA:** FAM-channel

***Alicyclobacillus acidocaldarius*-DNA:** CY5-channel

**Inhibition Control-DNA:** ROX-channel


A sample is **Alicyclobacillus spp** positive, if there is a detectable fluorescence increase in the **HEX-channel** and the negative controls show no amplification. The positive controls should have a positive fluorescence signal. The Inhibition Control in the ROX-channel may be positive or negative (depending on the amount of DNA or inhibitors in the sample reaction). In the negative controls it has to be positive.

A sample is ***Alicyclobacillus acidoterrestris*** positive, if there is a detectable fluorescence increase in the **FAM-channel** and the negative controls show no amplification. The positive controls should have a positive fluorescence signal. The Inhibition Control in the ROX-channel may be positive or negative (depending on the amount of DNA or inhibitors in the sample reaction). In the negative controls it has to be positive.

A sample is ***Alicyclobacillus acidocaldarius*** positive, if there is a detectable fluorescence increase in the **CY5-channel** and the negative controls show no amplification. The positive controls should have a positive fluorescence signal. The Inhibition Control in the ROX-channel may be positive or negative (depending on the amount of DNA or inhibitors in the sample reaction). In the negative controls it has to be positive.

A sample is negative, if there is no detectable fluorescence increase in the different channels and the positive controls have a positive fluorescence signal. The negative controls show no amplification. The Inhibition Control in the ROX-channel has to be positive in the sample and in the negative controls, a false negative result due to inhibitory effects is then excluded.

# analysis flowchart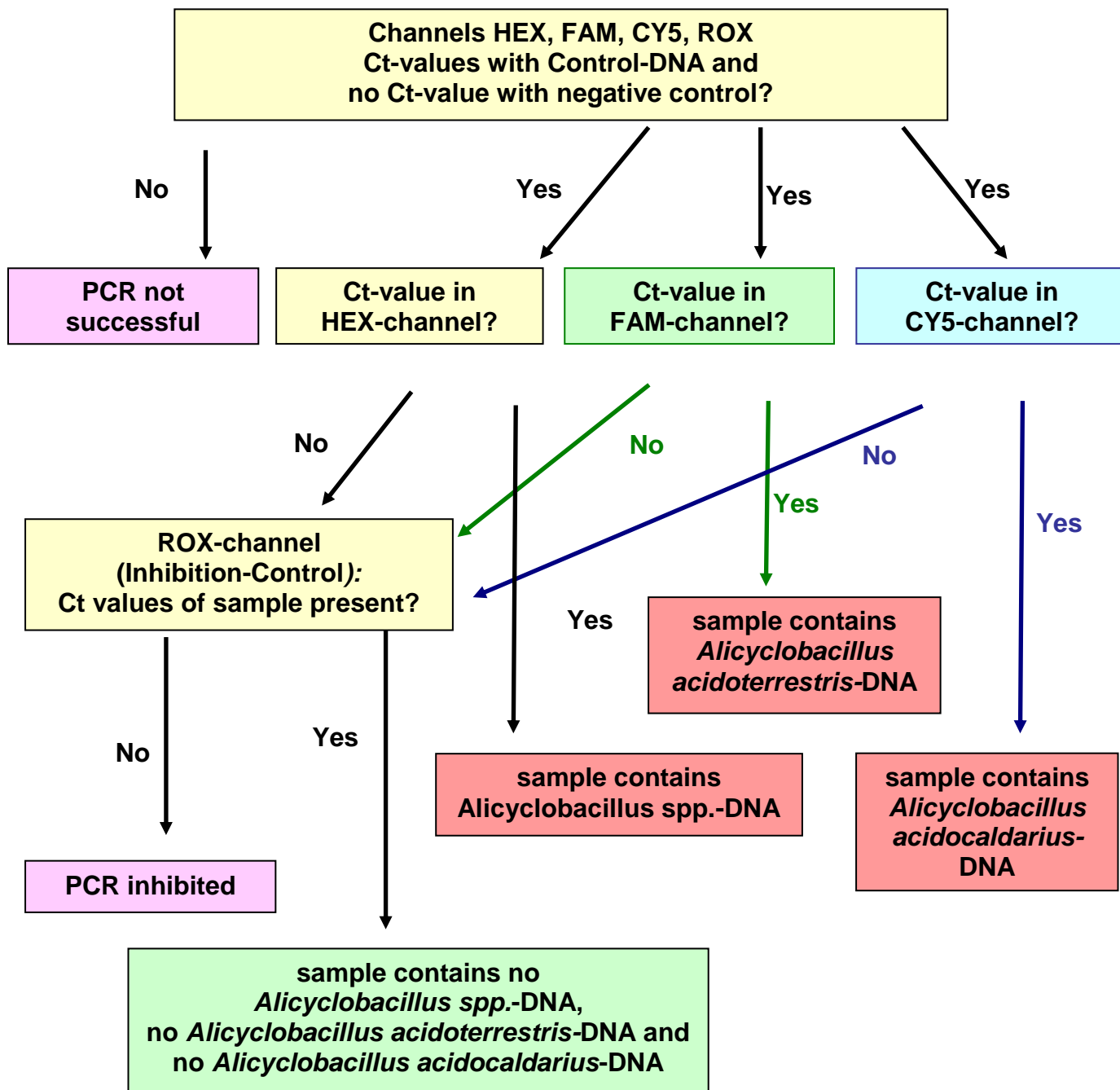

**Note:**

The polymerase-chain reaction (PCR) is protected by patents and requires a licence from Hoffmann-LaRoche Inc.. The provided product does not authorise the purchaser for the commercial use of this method.

GEN-IAL makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. If any materials are defective, GEN-IAL will provide a replacement product. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. GEN-IAL shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.