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Instructions For Use

Sample Buffer for bioavid allergen tests

Intended Use

The buffer is particularly suitable for the preparation of difficult samples (ketchup, mayonnaise) for the bioavid rapid allergen tests.

Storage

Store the buffer at room temperature. Do not use beyond the expiration date printed on the label.

Kit Contents

The Sample Buffer is available for 10 determinations, Art. No.: BS 815-10

815-100	Contents
1	Bottle with 100 ml Sample Buffer

Procedure

- Step 1: Add 9 ml Sample Buffer into a clean vial. Carefully close the bottle with remaining buffer.
- Step 2: Thoroughly ground and homogenize the sample.
- Step 3: Add 1 g or 1 ml sample into the vial with Sample Buffer. Close the vial.
- Step 4: Mix carefully avoiding foam formation (top down rolling) for 5 min.
- Step 5: Remove insoluble material (centrifuge for 1 min at 1000 x g).

Add 0.2 ml Running Buffer, then 0.1 ml of the prepared sample into a Reaction Vial of the selected bioavid allergen detection kit. Incubate the Reaction Vial with the sample for 5 min, tipping the vial a few times onto the work bench in order to mix antibody in the vial with the sample. After 5 min, insert the appropriate allergen test strip. Read the result as printed in the corresponding instruction leaflet.

Limitation

It is advised to run a spiked sample in parallel.

Examples for Detection Limits

Only approximate data can be given, as samples will differ significantly depending on composition and processing. Data are derived from spiking experiments.

Test kit	Allergen	Limit of Detection [ppm]		
		Mayonnaise	Ketchup	Fruit Juice
BL 613	Milk Powder	10	10	10
BL 607	Walnut	50		
BL 608	Egg yolk (uncooked)	25		

Note

Bioavid offers to their customers a chargeable service to analyze samples with respect to sample preparation and detection limit. Please contact your sales representative.