r-biopharm®

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# RIDA<sup>®</sup> Extraction Solution (colorless)

## 1. Intended use

### 1.1. Use with RIDASCREEN<sup>®</sup>FAST Casein (Art. No. R4612)

The RIDA<sup>®</sup> Extraction Solution (colorless) (Art. No. R7098) can be used for the sample preparation of sausages and bakery goods in the RIDASCREEN<sup>®</sup>FAST Casein test kit.

The sample preparation is described in the instruction for use or application of the test.

### 1.2. Use with Gliadin test systems

The RIDA<sup>®</sup> Extraction Solution (colorless) (Art. No. R7098) can be used for the sample preparation of raw materials, heated and processed food for the following qualitative dip stick tests:

- RIDA<sup>®</sup>QUICK Gliadin (Art. No. R7003)
- RIDA<sup>®</sup>QUICK Gliadin (single packaged) (Art. No. R7004)

It is recommended to use the Cocktail (patented) (Art. No. R7006 / R7016) for quantitative analysis with RIDASCREEN<sup>®</sup> Gliadin (Art. No. R7001) or RIDASCREEN<sup>®</sup>FAST Gliadin (Art. No. R7002).

The RIDA<sup>®</sup> Extraction Solution (colorless) should only be used after comparison with Cocktail (patented), because extraction with RIDA<sup>®</sup> Extraction Solution (colorless) may result in reduced recoveries in some samples. An application can be requested at R-Biopharm.

## 2. Reagent provided

1 x 105 ml Extraction Solution

## 3. Warnings and precautions for the users

The RIDA<sup>®</sup> Extraction Solution (colorless) should only be used by trained laboratory personnel. The instruction for use must be strictly followed.

The RIDA<sup>®</sup> Extraction Solution (colorless) is harmful to health. It contains mercaptoethanol. It should be worked under a chemical hood and skin contact should be avoided (use gloves).

Please refer to the component safety information in the material safety data sheet (SDS) for this product, available online at [www.r-biopharm.com](http://www.r-biopharm.com).

All reagents and materials must be recovered or disposed after use at customers own responsibility according to the protection of human health and the environment. Please observe the applicable national regulations concerning waste disposal (e.g. Waste Management Act, Regulations on Dangerous Chemicals, etc.).

#### 4. Storage instructions

Store the RIDA<sup>®</sup> Extraction Solution (colorless) at room temperature (20 - 25 °C / 68 - 77 °F). Please do not freeze.

Do not use the RIDA<sup>®</sup> Extraction Solution (colorless) after the expiration date (see test kit label).

**Further product information and applications, please contact your local distributor or R-Biopharm at this address: [sales@r-biopharm.de](mailto:sales@r-biopharm.de).**

#### Version overview

Version number	Chapter and title
2014-07-09	Release version
2021-09-13	General revision: <ul style="list-style-type: none"><li>– Adaptation of intended use in chapter 1</li><li>– Addition of proper disposal in chapter 3</li><li>– Additional sections "Version overview" and "Explanation of symbols"</li><li>– New disclaimer</li></ul>

## Explanation of symbols

General symbols:

	Follow the instruction for use
	Batch number
	Expiry date (YYYY-MM)
	Storage temperature
	Article number
	Number of test determinations
	Manufacturing date (YYYY-MM)
	Manufacturer + address

## Disclaimer

The user assumes all risk in using R-Biopharm AG's products and services.

R-Biopharm AG will warrant that its products and services meet all quality control standards set by R-Biopharm AG, and R-Biopharm AG will, at its option, replace or repair any components, product or repeat services which prove to be defective in workmanship or material within product specific warranty periods or expiration dates and which our examination shall disclose to our satisfaction to be defective as such.

This warranty is expressly in lieu of all other warranties, expressed or implied, as to quality, description, fitness for any particular purpose, merchantability, productiveness, or any other matter. R-Biopharm AG shall be in no way responsible for the proper use of its products and hereby disclaims all other remedies, warranties, guarantees or liabilities, expressed or implied, arising by law or otherwise, and it shall have no liability for any lost profits or damage, direct, indirect or otherwise, to person or property, in connection with the use of any of its products or services.

This warranty shall not be extended, altered or varied except by a written instrument signed by an authorized representative of R-Biopharm AG.

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are registered trademarks of R-Biopharm AG.