**CONGEN**

# **SureFood® QUANTARD Allergen 40**

Art. No. S3301

## **User Manual**

**April 2022**

 **Inhalt**

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## 1 General Information

### 1.1 Description

The SureFood® QUANTARD Allergen 40 is validated for quantitative determination of 1 mg to 400 mg allergenic substance / kg food sample.

The laboratory reference material SureFood® QUANTARD Allergen 40 is based on **corn flour** material. It contains all potentially allergenic food ingredients according to regulation (EU) No 1169/2011 (except for **molluscs, lactose and sulphur dioxide**) in a concentration of **40 mg/kg**.

#### Content of allergenic substances:

Allergen	Material
Gluten containing cereals	Wheat flour type 405 ( <i>Triticum</i> spp.)
Crustacean	Black Tiger shrimps ( <i>Penaeus monodon</i> )
Egg	Egg powder , hen ( <i>Gallus gallus</i> )
Fish	Humpback salmon ( <i>Oncorhynchus gorbuscha</i> )
Peanut	Peanut ( <i>Arachis hypogaea</i> ), untreated, peeled
Soy	Soybean ( <i>Glycine max</i> )
Milk	Milkpowder, cow ( <i>Bos taurus</i> )
Edible Nut	Hazelnut ( <i>Corylus avellana</i> ), peeled Almond ( <i>Prunus dulcis</i> ), untreated Walnut ( <i>Juglans regia</i> ), untreated Cashew ( <i>Anacardium occidentale</i> ), untreated Pecan ( <i>Carya illinoensis</i> ), untreated Brazil nut ( <i>Bertholletia excelsa</i> ), peeled Pistachio ( <i>Pistacia vera</i> ), peeled Macadamia ( <i>Macadamia ternifolia</i> ), roasted
Celery	Celery seeds ( <i>Apium graveolens</i> )
Mustard	Mustard seeds, yellow ( <i>Sinapis alba</i> )
Sesame	Sesam seeds ( <i>Sesamum indicum</i> )
Lupin	Lupin seeds ( <i>Lupinus hartwegii</i> ), peeled
Buckwheat	Buckwheat ( <i>Fagopyrum esculentum</i> )

The laboratory reference material SureFood® QUANTARD Allergen 40 is validated in combination with the DNA extraction method SureFood® PREP Advanced, protocol 1, and the SureFood® ALLERGEN real-time PCR detection methods.

Further information about the quantification of allergenic substances is given in the manual of the SureFood® ALLERGEN real-time PCR products.

In case of using alternative methods the sample weight and extraction method in combination with the detection system have to be validated.

**1.2 DNA-preparation**

For DNA-preparation the use of SureFood® PREP Advanced (Art. No. S1053), protocol 1 is recommended.

**1.3 Kit components and storage**

Reagent	Amount
QUANTARD Allergen 40	1 x 2.0 g

**Store all reagents dry at –20°C and protected from light.**

**Note: To avoid moisture absorption thaws the unopened laboratory reference material at room temperature for at least one hour before use.**

**1.4 Additionally required equipment and materials**

- DNA-Extraction kit (e.g. SureFood® PREP Advanced, Art. No.: S1053)
- Real-time PCR methods (SureFood® ALLERGEN series)

**2 Further Information****2.1 Technical Support**

For further questions please contact your distributor or send an e-mail to [sales@r-biopharm.de](mailto:sales@r-biopharm.de).

**2.2 Distribution and Ordering**

R-Biopharm AG  
An der neuen Bergstrasse 17,  
64297 Darmstadt, Germany  
Phone: +49 (0) 61 51 - 81 02-0  
Fax: +49 (0) 61 51 - 81 02-20  
E-Mail: [orders@r-biopharm.de](mailto:orders@r-biopharm.de)  
[www.r-biopharm.com](http://www.r-biopharm.com)

r-biopharm®