**CONGEN**

# **SureFood® ALLERGEN Pine Nut**

Art. No. S3624  
100 rxn

## **User Manual**

**January 2022**

 **Inhalt**

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## **Content**

|       |   |    |
|-------|---|----|
| 1     | General Information .....                           | 9  |
| 1.1   | Description .....                                   | 9  |
| 1.2   | Limit of Detection .....                            | 9  |
| 1.3   | DNA-preparation .....                               | 10 |
| 1.4   | Kit components and storage .....                    | 10 |
| 1.5   | Additionally required equipment and materials ..... | 10 |
| 1.6   | Setup .....   | 11 |
| 1.7   | Detection channel Set-up .....                      | 11 |
| 2     | Qualitative Analysis .....                          | 12 |
| 2.1   | Protocol .....                                      | 12 |
| 2.1.1 | Preparation of the master-mix .....                 | 12 |
| 2.1.2 | Preparation of the real-time PCR-mix .....          | 12 |
| 2.2   | Interpretation of results .....                     | 13 |
| 3     | Further Information .....                           | 14 |
| 3.1   | Product Information .....                           | 14 |
| 3.2   | Technical Support .....                             | 14 |
| 3.3   | Distribution and Ordering .....                     | 14 |

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## **1 General Information**

### **1.1 Description**

The SureFood® ALLERGEN Pine Nut is a real-time PCR for the direct, qualitative detection of specific pine nut (*Pinus* spp.) DNA sequences in food.

Each reaction contains an internal amplification control (IAC). If the DNA contains PCR inhibiting substances, the signal of the amplification control will be affected or the amplification will be suppressed. Examples for PCR inhibiting substances are alcohols (e.g. ethanol, isopropanol), surfactants (e.g. CTAB, SDS, Triton X100) and salts (e.g. sodium chloride). In addition spices, herbs, algae, cocoa and further sample matrices might have PCR inhibiting effects.

The real-time PCR assay can be performed with commonly used real-time PCR instruments, equipped for detection of two fluorescence emissions at the channels FAM and VIC/HEX at the same time. The technical validation of instruments was performed on Agilent AriaDx, Bio-Rad CFX96 Dx, R-Biopharm RIDA®CYCLER, Roche LightCycler® 480 II and Qiagen Rotor-Gene Q.

### **1.2 Limit of Detection**

The SureFood® ALLERGEN Pine Nut PCR has a limit of detection of  $\leq 1$  mg / kg using SureFood® PREP Advanced, protocol 1.

The limit of detection of the complete method (DNA extraction and real-time PCR) depends on sample matrix, processing grade, DNA-preparation and DNA-content.

The SureFood® PCR systems are very sensitive and therefore even a small amount of target DNA is sufficient for a successful analysis. The concentration of total DNA in the sample does not allow a conclusion on the quantity and quality of the target DNA.

**1.3 DNA-preparation**

For DNA-preparation the use of SureFood® PREP Advanced (Art. No. S1053), protocol 1 or SureFast® Mag PREP Food (Art. No. F1060) is recommended.

**1.4 Kit components and storage**

| <b>Kit Code</b> | <b>Reagent</b>   | <b>Amount</b> | <b>Lid Color</b> |
|-----------------|------------------|---------------|------------------|
| 1               | Reaction Mix     | 2 x 1100 µl   | Yellow           |
| 2               | Taq Polymerase   | 1 x 80 µl     | Dark Red         |
| 3               | Positive Control | 1 x 200 µl    | Light Blue       |

**Store all reagents at –20°C and protected from light. The Taq Polymerase can be stored at +2 to +8°C for multiple uses on the same day.**

**Note: The Taq Polymerase may be in a frozen or unfrozen state. This does not affect the quality of the Taq Polymerase or the performance of the real-time PCR.**

**1.5 Additionally required equipment and materials**

- DNA-Extraction kit (e.g. SureFood® PREP Advanced Art. No. S1053 or SureFast® Mag PREP Food Art. No. F1060)
- real-time PCR instrument with two detection channels (510 nm and 580 nm)
- real-time PCR consumables (plates, tubes, foils, caps)
- pipettes with filter tips
- powder-free disposable gloves
- Vortex mixer
- micro centrifuge with a rotor for the reaction tubes

**1.6 Setup**

|  | <b>Blockcyler &amp;<br/>R-Biopharm RIDA®CYCLER</b> | <b>Rotorcyler</b> |
|--|--|-------------------|
| Initial Denaturation (HOLD)            | 5 min, 95°C  | 1 min, 95°C       |
| Cycles                                 | 35   | 35                |
| Denaturation                           | 15 sec, 95°C                                       | 10 sec, 95°C      |
| Annealing/Extension (CYCLE)            | 30 sec, 60°C                                       | 15 sec, 60°C      |
| Temperature Transition Rate/ Ramp Rate | Maximum  | Maximum           |

**1.7 Detection channel Set-up**

| <b>Real-time PCR device</b>              | <b>Detection</b> | <b>Detection channel</b> | <b>Quencher</b> | <b>Note</b>  |
|--|------------------|--------------------------|-----------------|--|
| <b>Agilent<br/>AriaDx/Mx</b>             | pine nut         | FAM                      | +               |  |
|  | IAC              | HEX                      | +               |  |
| <b>Bio-Rad<br/>CFX96/Dx</b>              | pine nut         | FAM                      | +               |  |
|  | IAC              | VIC/HEX                  | +               |  |
| <b>R-Biopharm<br/>RIDA®CYCLER</b>        | pine nut         | green                    | +               |  |
|  | IAC              | yellow                   | +               |  |
| <b>Qiagen Rotor-<br/>Gene Q</b>          | pine nut         | green                    | +               | <b>Note:</b> Please use only 0.1 ml reaction tubes. The gain settings must be set to 5 (factory default) for all channels. |
|  | IAC              | yellow                   | +               |  |
| <b>Roche<br/>LightCycler®<br/>480 II</b> | pine nut         | 465-510                  | +               |  |
|  | IAC              | 533-580                  | +               |  |

## 2 Qualitative Analysis

### 2.1 Protocol

#### 2.1.1 Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions) for the specific PCR assay.

Recommended control reactions for the specific PCR assay: negative control, extraction control, positive control. The reaction mix contains an internal amplification control (IAC) per reaction.

#### Reactions needed for the qualitative pine nut detection:

3 reactions for controls (1x no-template control, 1x extraction control, 1 positive control)

For each sample: at least 1 reaction for each sample DNA

It is also recommended to prepare the master-mix with 10 % additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix and centrifuge before opening and use.

#### Example for the calculation and preparation of 10 reactions:

| Components of the master-mix | Amount per reaction | 10 reactions (with 10% excess) |
|------------------------------|---------------------|--------------------------------|
| Reaction Mix                 | 19.3 µl             | 212.3 µl                       |
| Taq Polymerase               | 0.7 µl              | 7.7 µl                         |
| <b>Total volume</b>          | <b>20 µl</b>        | <b>220 µl</b>                  |

**Mix each master-mix well and centrifuge shortly before use.**

#### 2.1.2 Preparation of the real-time PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells.
- Close the negative control (the negative control is ready for PCR without any addition).
- Pipette 5 µl of sample DNA into the designated tubes/wells and close them.
- Pipette 5 µl of Positive Control into the designated tubes/wells and close them.
- Centrifuge all tubes/plates shortly at low speed.
- Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

## **2.2 Interpretation of results**

The evaluation has to be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer.

The control reactions have to show the correct results.

Pine nut DNA is detected in the FAM-channel. In the VIC/HEX-channel the amplification control is detected.

A sample is stated **positive**, if the sample DNA shows amplification in the detection system. High amplicon concentrations can result in a weak or absent signal of the internal amplification control (IAC).

A Cp value for the internal amplification control (IAC) is not needed to obtain a positive result of the positive control.

A sample is stated **negative**, if the sample DNA shows no amplification in the detection system and if the internal control (VIC/HEX-channel) of the sample is **positive** with a shift in Cp-value  $\leq 2$  compared to the negative control.

If the sample DNA in the VIC/HEX-Channel shows **no amplification** or a shift in Cp-value  $> 2$  compared to the negative control, it contains PCR inhibiting substances. A significant decrease in the fluorescence signal can also show the presence of PCR inhibiting substances. Under these circumstances DNA isolation and purification of the sample need to be improved. Alternatively the DNA can be diluted (recommendation 1:2 in PCR-water) and analysed again for inhibition. Please note that the dilution factor also affects the detection limit of the specific pine nut PCR assay.

Alternatively, a self-selected reference sample at the detection limit can be used to assess the sample. This reference sample is used outside CONGEN's quality and evaluation guidelines.

### **In general:**

Samples tested negative still could contain an allergen contamination below the limit of detection of the assay, or they might contain other allergen components like proteins or lipids for example.

### **3 Further Information**

#### **3.1 Product Information**

- Detailed information about setup of several real-time PCR devices  
(Download: [www.congen.de/en/company/downloads](http://www.congen.de/en/company/downloads))
- Validation Report upon request

#### **3.2 Technical Support**

For further questions please contact your distributor or send an e-mail to [sales@r-biopharm.de](mailto:sales@r-biopharm.de).

#### **3.3 Distribution and Ordering**

R-Biopharm AG  
An der neuen Bergstrasse 17,  
64297 Darmstadt, Germany  
Phone: +49 (0) 61 51 - 81 02-0  
Fax: +49 (0) 61 51 - 81 02-20  
E-Mail: [orders@r-biopharm.de](mailto:orders@r-biopharm.de)  
[www.r-biopharm.com](http://www.r-biopharm.com)